

# CAPP Origami

## REAGENT RESERVOIRS



## Fold, pour and pipette...

The **CAPP Origami™** represents an innovative way of thinking when it comes to reagent reservoirs.

### Perfect V-shape maximises the full retrieval of all reagents

The shape of the reservoir, along with the stiffness of the material used, ensures that the reservoir forms a sharp V-shape at the bottom. This makes your pipetting easier and also enables you to save on reagents as you only have to pour the amount needed into the reservoir.

### Flick-up design saves space when stored

**CAPP Origami™** reservoirs are packed as flat sheets which means they take up considerably less room in storage.

### Easy to use

To use a **CAPP Origami™** reservoir, simply flick it up by pushing the sides together. This is very easy due to the pre-folded lines. As soon as the **CAPP Origami™** reservoir reaches the proper shape you can pour liquid in and start pipetting straight away. The folds and the polypropylene material ensure that the reservoirs stay upright once folded into the correct reservoir shape.


## We take pipetting to perfection.



Take flat reservoir from pack



push two outer sides towards the middle

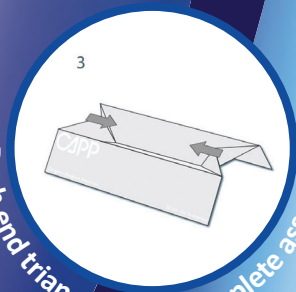

push end triangles in to complete assembly

# Ordering Information

Cat. No.	Description
CA40505	CAPPOrigami 30mL (8- and 16-channel pipettes), bag w/ 50 pcs.
CA40506	CAPPOrigami 30mL (8- and 16-channel pipettes), Pre-sterile, 10 bags w/ 5 pcs. each
CA40510	CAPPOrigami 40mL (12-channel pipettes), bag w/ 50 pcs.
CA40511	CAPPOrigami 40mL (12-channel pipettes), Pre-sterile, 10 bags w/ 5 pcs. each

## CAPP*Aero 96* Multichannel Pipettes

Cat. No.	Description	Inaccuracy %	Imprecision %	Colour
C10-8	CAPPAero 96 multi pipettes, 8-channel, 0.5-10µL	5.00 /2.00	3.60 /1.00	○
C20-8	CAPPAero 96 multi pipettes, 8-channel, 2-20µL	5.00 /2.00	3.60 /1.00	●
C50-8	CAPPAero 96 multi pipettes, 8-channel, 5-50µL	3.20 /1.40	1.60 /0.60	●
C100-8	CAPPAero 96 multi pipettes, 8-channel, 10-100µL	2.50 /1.30	1.20 /0.50	●
C200-8	CAPPAero 96 multi pipettes, 8-channel, 20-200µL	1.80 /1.20	0.80 /0.40	●
C300-8	CAPPAero 96 multi pipettes, 8-channel, 30-300µL	1.66 /1.40	0.60 /0.30	●
C10-12	CAPPAero 96 multi pipettes, 12-channel, 0.5-10µL	5.00 /2.00	3.60 /1.00	○
C20-12	CAPPAero 96 multi pipettes, 12-channel, 2-20µL	5.00 /2.00	3.60 /1.00	●
C50-12	CAPPAero 96 multi pipettes, 12-channel, 5-50µL	3.20 /1.40	1.60 /0.60	●
C100-12	CAPPAero 96 multi pipettes, 12-channel, 10-100µL	2.50 /1.30	1.20 /0.50	●
C200-12	CAPPAero 96 multi pipettes, 12-channel, 20-200µL	1.80 /1.20	0.80 /0.40	●
C300-12	CAPPAero 96 multi pipettes, 12-channel, 30-300µL	1.66 /1.40	0.60 /0.30	●



Uthleber Weg 14  
99734 Nordhausen  
Germany  
Tel. +49 (0) 3631 65242-0  
Fax. +49 (0) 3631 65242-90  
info@ahn-bio.com  
[www.ahn-bio.com](http://www.ahn-bio.com)  
V. 2024-12

Distributed by:

